

GRS Hotstart Taq DNA Polymerase
#GE71.0500 | # GE71.2500 | # GE72.0500 | # GE72.2500
 (for research only)

- Product:** GRS Hotstart Taq DNA polymerase consists of GRS Taq bound to a specific antibody that blocks polymerase activity, preventing or minimizing the synthesis of non-specific products and primer-dimers prior to PCR cycling, which results in an overall higher specificity. The heat-labile antibody can be rapidly inactivated by increasing the temperature for a few minutes to 95°C (hot-start) resulting in polymerase activation. Like GRS Taq, this enzyme has 5'→3' polymerase activity but no 3'→5' exonuclease activity (no proofreading). It is suitable for amplifying DNA targets up to 5kb and produces 3' A-overhangs allowing TA-cloning
- Source:** Purified from *E.coli* harbouring the DNA polymerase gene from *Thermus aquaticus* YT-1.
- Quantity:** **#GE71.0500** and **#GE72.0500** comprises 1 vial of 500U of GRS Hotstart Taq, **#GE71.2500** and **#GE72.2500** comprises 5 vials of 500U of GRS Hotstart Taq, at a concentration of 5U/μl in 20mM Tris-HCl (pH 8.0), containing 0.1mM EDTA, 10mM β-mercaptoethanol, 222mM ammonium sulphate, 0.5% (v/v) Triton X-100, 0.5% (v/v) Tween®20, and 50% glycerol.
- Contents:** GRS Hotstart Taq Polymerase is supplied with PCR Buffer (10x concentrated), consisting of 100mM Tris (pH 8.9), 500mM KCl, supplemented with BSA (2.5mg/ml) and with a separate vial of 25mM MgCl₂.

#GE72.0500 and **#GE72.2500** are also supplied with separated dNTPs.

Product	GE71.0500	GE71.2500	GE72.0500	GE72.2500
GRS Hotstart Taq (5U/μl)	100μl	5x 100μl	100μl	5x 100μl
PCR Buffer (10x)	1,5ml	5x 1,5ml	1,5ml	5x 1,5ml
MgCl ₂ (25mM)	1,5ml	5x 1,5ml	1,5ml	5x 1,5ml
dNTPs (10mM each)	-	-	1,0ml	3x 1,0ml

- Applications:** Routine PCR
- Properties:** Amplicon size: up to 5kb.
 Extension Rate 1kb/min
 Hotstart: Yes
 A-overhang: Yes
- QC:** Functionally tested in PCR. Bacterial DNA free (tested by PCR). Absence of endonucleases, exonucleases, and ribonucleases was confirmed by appropriate assays.
- Storage:** -20°C for at least 1 year.

GRiSP Research Solutions
 Rua Alfredo Allen, 455
 4200-135 Porto
 Portugal
www.grisp.pt | info@grisp.pt

Basic Protocol

Optimal PCR cycling conditions (incubation times, temperatures, and concentrations) depend on DNA target (GC-content, size, quantity, purity, etc), specific primers, and buffer composition, MgCl₂-concentration and GRS Hotstart Taq-concentration, and need to be determined case by case. We suggest to start with the following basic protocol and subsequently optimize MgCl₂-concentration, annealing temperature, incubation times and cycling number.


Mix for each PCR reaction, starting with the greatest volume (usually water) and ending with GRS HotstartTaq:

Component	Volume	Final Conc.
MgCl ₂ (25mM)	2,5 µl	2,5 mM
dNTPs (10 mM each)	1 µl	0,4 mM
Forward primer (5 pmol/ µl)	1,5 µl	0,3 µM
Reverse primer (5 pmol/ µl)	1,5 µl	0,3 µM
PCR Buffer (10x)	2,5 µl	1X
Template DNA	0,25 - 10 µl	2-200 ng/µl*
GRS Hotstart Taq (5U/ul)	0,25 µl	0.05U/µl
PCR -grade water	up to 25 µl	

*) In case of plasmid DNA use 1-3 ng, in case of gDNA add 200-300ng
For smaller/larger reaction volumes, scale it down/up proportionally.

In order to minimize risk of contamination, reagent loss and improve pipetting accuracy, we recommend to prepare a mastermix for multiple samples (N), always including a negative control for the detection of possible contaminants, by mixing all components (N+1), except template DNA, dividing the mixture equally into each tube and then add template DNA or in PCR grade water in case of the control to the individual PCR tubes.

Set-up PCR amplification as follows:

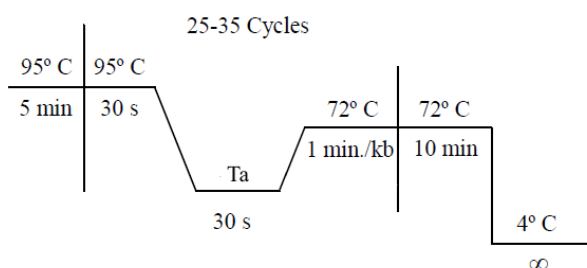

After an initial cycle of 5 min at 95°C (antibody inactivation and complete denaturation of template DNA including removal of all secondary DNA structures such as hairpins), cycle 25-35 times for 30 seconds at 95°C, 30 seconds at Ta and 1 min (per kb of target DNA) at 72°C for extension. Set the annealing temperature (Ta) as the melting temperature (Tm) of the primer with the lowest Tm. After amplification, include a final extension step of 10 min at 72°C to ensure that all amplicons are fully extended and include 3'-A-overhang. Analyze PCR products by DNA Agarose gel electrophoresis.

Optimization

Annealing Temperature (Ta)

Optimizing the annealing temperature is crucial, as a too low temperature might result in non-specific amplification whereas a too high temperature results in no amplification. The melting temperature (Tm) is defined as the temperature in which 50% of the primer and its complementary sequence of the target DNA are present as duplex DNA. By increasing the temperature above the melting temperature, this percentage decreases, however, primers will still ligate (up to a certain point) and initiate extension. PCR can therefore be performed at temperatures of several degrees higher than Tm and it is therefore recommended to optimize the Ta by performing a temperature gradient.

MgCl₂

GRS Hotstart Taq DNA Polymerase is supplied with a PCR Buffer (10x) without magnesium and a separated vial of 25mM MgCl₂. Taq requires MgCl₂, however, the final concentration may have great influence on both quantity and specificity of the amplification and we highly recommend to optimize the final MgCl₂-concentration (normally between 1,5 and 3,5mM in steps of 0.5mM).

Incubation times

Denaturation and annealing steps may require less time depending on the thermocycler apparatus, reaction volume and PCR tube (varies with the efficiency of heat-transfer). It might be worthwhile to optimize (reduce) times to as low as 10-20 seconds for both denaturation and annealing steps, which will greatly reduce overall PCR time.

GC-rich targets

If having troubles amplifying GC-rich targets, one could try adding 5-10% DMSO, up to 10% glycerol, or 1-2% formamide (or any combination of these solvents) to the reaction mixture. Note that the optimal Ta will decrease approximately 5°C in 10% DMSO.